



## Emergency Shower Station Weekly Inspection Checklist

**How to Test:**

1. Is the location of the emergency equipment well lit and identified with a highly visible sign?
2. Is access to the shower free of obstructions from all areas of the laboratory or hallway?
3. Flush shower for at least 10 seconds. Place a 5-gallon bucket underneath the safety shower and a shower curtain around the shower head and inside the bucket to catch all the water.
4. Is the water clear?
5. Evaluate for adequate flow.
6. Does the water continue to flow until the pull handle is turned off?
7. Is the plumbing free from leakage?
8. Promptly clean up any water that is spilled on the floor as a result of this process. Any water collected can be emptied at the most convenient location such as a sink, custodial floor sink or a floor drain.
9. Date and initial the weekly inspection log below.

**How to obtain assistance:**

**Questions, maintenance requests, training information:**

EH&S 575-5448

**Maintenance requests may also be called in directly:**

Service Desk: 575-5050

Facilities Services (Zone Maintenance):

Zone A: 575-3730 Zone B: 575-6173

Zone D: 575-3414 Zone E: 575-2968

Zone F: 575-7354

**Emergencies:**

Immediately stand in the water stream and flush the affected area for 15 minutes.

Remove contaminated clothing while showering and not before. CALL 911 and seek medical assistance.

Date	Initials	Comment		Date	Initials	Comment		Date	Initials	Comment